	记录编号 FN: SMP-RD-A0002-R06	版本号 Ver: 2.1
	归档 Filing: filed immediately, permanently kept	生效日期 Ed: 2023.01.12

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多宁/DuoNing

动物细胞高性能培养基 High-Performance Culture Medium for Animal-Cells

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The high-efficiency feed medium B is used together with 2×103A high-efficiency feed medium A, 2×104 high-efficiency feed medium A, 2×105 high-efficiency feed medium A, 2×107 high-efficiency feed medium A and 103 high-efficiency feed medium A. Among them, 2×103A high-efficiency feed medium is a kind of high-efficiency fed-batch culture medium without animal-derived components and protein, which is suitable for the research and development of therapeutic protein products by Chinese hamster ovary (CHO) and the fed-batch culture in the production process. This product contains no hypoxanthine, thymine, L- glutamine, P188 and sodium bicarbonate. It is suitable for the culture of different cell lines, such as CHO-K1, CHO-S, DG44 and CHOZN, using GS system and DHFR screening system. 2×103A high-efficiency feed medium is divided into 2×103A high-efficiency feed medium A and high-efficiency feed medium B.

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2×103A high-efficiency feed medium is an efficient feeding formula specially designed to improve the growth and production performance of CHO cells. It is rich in nutrients such as amino acids and vitamins, which can meet the nutritional needs of high-density culture of cells and high expression of products. Combined with the characteristics of cell

	记录编号 FN: SMP-RD-A0002-R06	版本号 Ver: 2.1
	归档 Filing: filed immediately, permanently kept	生效日期 Ed: 2023.01.12

lines, reasonable feeding strategy can obviously improve the growth of cells and the antibody expression.

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It is suggested that the first feeding should be started in the middle and late period of cell exponential growth, when the cell density is above 2.0×10^6 cells/mL. The recommended feeding amount is: 5% on the 3rd day, 5% on the 5th day, 7.5% on the 7th day, 7.5% on the 9th day, 5% on the 11th day and 5% on the 13th day.

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It is suggested that the total feeding amount should not exceed 40%.

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$2 \times 103A$ high-efficiency feed medium can be used in combination with BR7, and the culture efficiency can be further improved. The recommended concentration of BR7 for each addition is 1%.

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Hypoxanthine and thymine were added as needed when cultivating GS screening system engineering cell lines.

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3~8mM L- glutamine was added as needed when cultivating DHFR screening system engineering cell lines.

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The liquid packaging of this product will partially precipitate when it is placed at 2~8°C, so it is recommended to use it within two months.

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Suitable for powder packaging (taking 1L as an example)

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Prepare ultrapure water (35 ~ 40°C) with 80% of the liquid volume;

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Add 99.03g of $2 \times 103A$ high-efficiency feed medium A powder, stir for 20 min, and dissolve completely. At this time, the solution is slightly turbid, and the pH is about 4.0;

- 另外准备 50ml 0.4M NaOH 溶液，加入高效补料培养基 B 粉末 2.96g，搅拌 10 min，溶解完全，记为 B 溶液；
In addition, prepare 50 ml of 0.4M NaOH solution, add 2.96g of high-efficiency feed medium B powder, stir for 10 min, and record it as solution B;
- 加入 B 溶液，继续搅拌 20 min；
Add solution B and continue stirring for 20 min;
- 调节 pH 至 6.70 左右，溶液澄清；
Adjust the pH to about 6.70, and the solution will be clarified;
- 加入 2.1g 碳酸氢钠，调节 pH 至 6.90~7.10；
Add 2.1g sodium bicarbonate to adjust the pH to 6.90 ~ 7.10;
- 定容，搅拌 5~10 min；
Constant volume, stirring for 5 ~ 10 min;
- 用 0.22 μ m 过滤器除菌过滤；
Sterilize and filter with a 0.22 μ m filter;
- 2 \times 103A 培养基液体渗透压范围在 750~850 mOsm/kg。
The osmotic pressure of 2 \times 103A medium is 750~850 mOsm/kg.

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Shanghai production base, powder packaging: 2 $^{\circ}$ C to 8 $^{\circ}$ C, protect from light; validity period: 24 months.

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Wuxi production base, powder packaging: 2 $^{\circ}$ C to 8 $^{\circ}$ C, protect from light; retest date: 24 months.

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地址 Address: 上海市松江区民强路 1525 弄 30 幢 Building 30, Lane 1525, Minqiang Road, Songjiang District, Shanghai

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